


PinAAcle Burner System Components

- 1** Burner Head
- 2** Burner Head O-Ring
- 3** Mixing Chamber
- 4** Flow Spoiler
- 5** End Cap Assembly
- 6** Stainless Steel Nebulizer
- 7** Retaining Ring
- 8** Burner Head Adjustment Tool

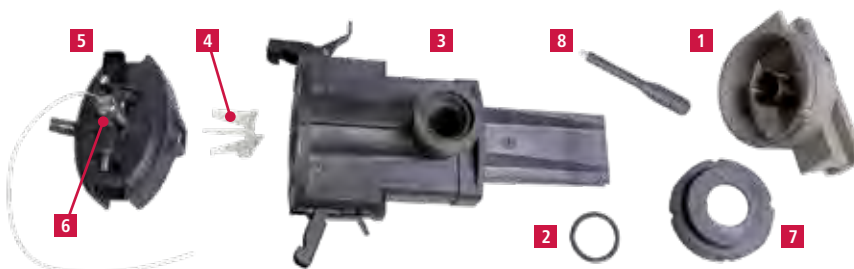

PinAAcle Burner System Components


Burner System Components

Description	PinAAcle 500/900 Part No.	AAAnalyst 100/200/300/400/700/800 Part No.
Viton Burner Head O-Ring Kit, 0.984 in. x 0.139 in. 09200079 O-Ring (3), 09200078 Super O-Lube (1)	09902219	09902219
Burner Mixing Chamber Assembly (Note: PinAAcle assembly includes components 2, 3, and 7; AAAnalyst component 2 only.)	N3160506	00570948
Burner Chamber Holder	N3160809	
Clamping Ring		N0401171
Flow Spoiler	00572561	00572561
Viton O-Ring for End Cap Gas Lines (Neb., Aux., Fuel) 0.312 in. x 0.070 in. (Qty. 4 required)	09200253	
Viton O-Ring for End Cap 1.892 in. x 0.103 in.	09902147	09902147
Kalrez O-Ring for End Cap Gas Lines (Neb., Aux., Fuel) 0.312 in. x 0.070 in. (Qty. 4 required)	09200275	
Kalrez O-Ring for End Cap 1.892 in. x 0.103 in.	09200277	09200277
Organics (Corkprene) O-Ring for End Cap	00472014	00472014
End Cap Assembly with Viton O-Rings (for use with metal body nebulizers)	N3160505	00570984
End Cap Assembly with Kalrez O-Rings (for use with metal body nebulizers)	N3160188	
End Cap Assembly with Viton O-Rings (for use with plastic body nebulizers)	N3160511	N0370392
End Cap Assembly with Kalrez O-Rings (for use with plastic body nebulizers)	N3160189	
End Cap Assembly Impact Bead (for use with metal body nebulizers)	N3160545	
Burner Gasket Kit for End Cap Assembly 00570984 ; 00472014 (5), 09902147 (3), 09902219 (1), 09902236 (3), 09902240 (3)		00470988
Drain Tubing – Tygon Tubing with FEP lining	02507987	02507987
Pyrex Impact Bead Used with End Caps for Metal Body Nebulizers	00572615	00572615
Seal Plug for the 00572615 Pyrex Impact Bead used with end caps for metal body nebulizers	N3161224	00572624
O-Ring for seal plug in end caps with Impact Bead (Qty. 2 required)	09902236	09902240
Burner Drain Assembly for Organic Solutions	N3160683 ***	N3150230 *
Burner Drain Assembly (includes tubing and drain interlock, drain bottle 09200486 not included)		N2011074 *
Burner Drain Assembly (includes tubing and drain interlock, drain bottle 09200486 not included)	N0400058 ***	N0400058 **
Wide Mouth 15 L Drain Bottle	09200486	09200486

*For AAAnalyst 700/800, **For AAAnalyst 200/400, ***Also compatible with AAAnalyst 100/200/300/400

AAAnalyst Burner System Components

- 1** Burner Head
- 2** Burner Mixing Chamber
- 3** Clamping Ring
- 4** Flow Spoiler
- 5** End Cap O-Ring
- 6** End Cap Assembly

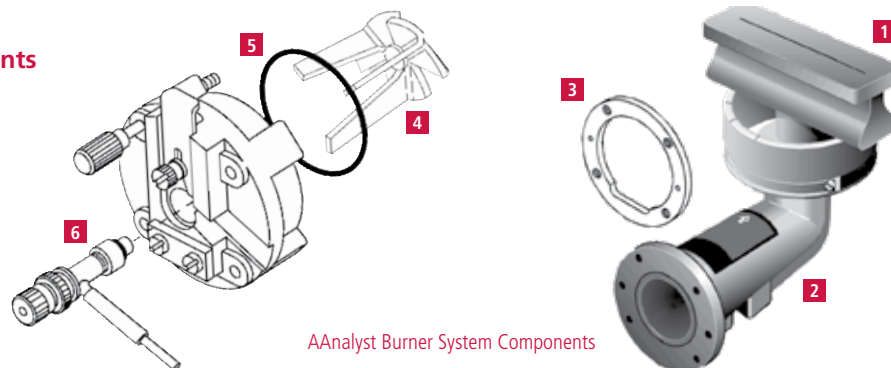

AAAnalyst Burner System Components