


Avio Injector Assembly

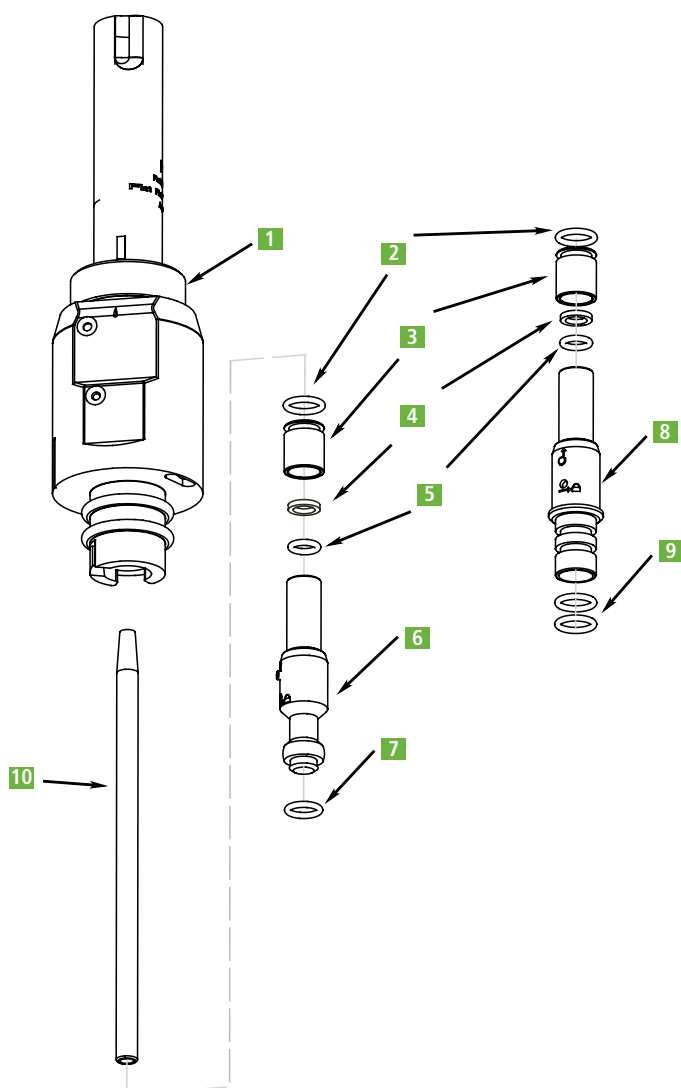

Description	Part No.
1 Knurled Nut	
2 O-Ring – Color Coded Blue	N0791333
3 Lock Collar	N0791163
4 Washer	N0776014
5 O-Ring	09902207
6 Cyclonic Injector Adapter	N0790143
7 O-Ring	09926070
8 Scott Injector Adapter	N0790154
9 O-Ring	09200518
10 Injector 2 mm Alumina	N0791183



Avio Injector Assembly Replacement Parts

Description	Part No.
Cyclonic Injector Assembly	N0790141
Scott Injector Assembly	N0790156
O-Ring Kit for Avio Scott Injector	N0790440
O-Ring Kit for Avio Cyclonic Injector	N0790441
O-Ring Kit for Avio Torch	N0790437
O-Ring Removal Tool	09220410

Injector Support Adapters and Replacement Parts

Injector Support Adapters

ICP Model	Spray Chamber	Part No.
Avio 200/500 Includes one of each: 09902207 , 09926070 , N0791333 , N0776014 , N0790143 , N0791163	Cyclonic	N0790141
Avio 200/500 Includes one of each: 09902207 , 09200518 , N0791333 , N0776014 , N0790154 , N0791163	Scott	N0790156
2x00/4x00/5x00/7x00 DV 8x00	All	N0770603
2x00/4x00/5x00/7x00 DV 8x00*	All	N0776091
4300V/5300V/7300V (bottom)	Cyclonic	N0771526
4300V/5300V/7300V (top)	Cyclonic	N0771527

*For 1.2 mm entire length injectors **N0776093** and **N0781007**

Injector Support Adapter O-Ring Kits

ICP Model	Part No.
Avio 200/500 Cyclonic	N0790441
Avio 200/500 Scott	N0790440
2x00/4x00/5x00/7x00 DV/8x00	N0770438
3000/3000 SCR/3X00 RL	N0690269
3x00 XL/3x00 DV/3000 SCX	N0690769