


TurboMatrix 40 Headspace Trap

Sample Head Assembly Replacement Parts

| Description | Part No. |
|---|-----------------|
| Platinum/Iridium Needle, Wide-bore | B0144169 |
| Platinum/Iridium Needle, Small-bore | B0500959 |
| Platinum/Iridium Needle, Jet | B0510364 |
| 1 Silcosteel Needle, for Headspace Trap Only | N6700130 |
| Stainless Steel Needle, Wide-bore | B0131385 |
| Stainless Steel Needle, Small-bore | B0500987 |
| Stainless Steel Needle, Jet (Ships with Instrument) | B4000011 |
| 2 Needle Seal Assembly (Without O-Rings) | B0500833 |
| 3 O-Ring for Needle Seal Assembly (pkg. 10) | B0198110 |
| 4 Vespel Ferrule 1/16 in. (pkg. 10) | 09920127 |
| 5 Male Nut 1/16 in. | N9302832 |
| 6 GLT Adapter Tube | B0503956 |
| GLT Adapter Tube, Silcosteel | N6700113 |
| 7 Graphite/Vespel Ferrule 1/16 in. x 0.4 mm For use with 0.25 mm ID Transfer Line (pkg. 10) | 09920104 |
| Graphite/Vespel Ferrule 1/16 in. x 0.5 mm For use with 0.32 mm ID Transfer Line (pkg. 10) | 09920105 |
| 8 Nut 1/16 in. Swagelok | N9300059 |
| 9 Fused-Silica Capillary Transfer Line: 0.25 mm ID x 5 m Length | N9301356 |
| 0.32 mm ID x 5 m Length | N9301357 |

Headspace Sample Head Assembly Schematic

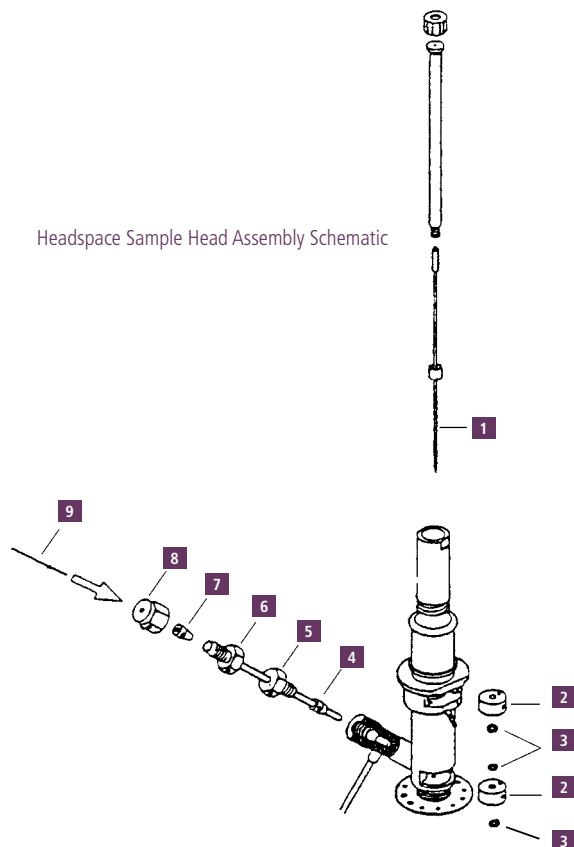

Solid Glass Blocking Trap

| Description | Part No. |
|--|-----------------|
| Block for Use in Standard Headspace Mode | N6701170 |

Sample Trays

For use on the Mid-Range or High-Capacity headspace sampler.

| Description | Part No. |
|---|-----------------|
| TurboMatrix 40 Mid-Range Sample Tray | M0413592 |
| TurboMatrix 110 High-Capacity Sample Tray | M0413593 |

Transfer Lines

| Description | Tubing ID (mm) | Length | Part No. |
|---------------------------------|----------------|--------|-----------------|
| Siltek Deactivated Fused Silica | 0.25 | 5 m | N9316607 |
| Siltek Deactivated Fused Silica | 0.32 | 5 m | N9316608 |

Miscellaneous Accessories

| Description | Part No. |
|---|-----------------|
| Gas Chromatography – Theory and Practice, Static Headspace Book by L. Ettre and B. Kolb | N1011210 |

Cold Trap Options

Headspace Trap instruments only.

| Description | Part No. |
|--|-----------------|
| TurboMatrix HS Trap Cold Trap Tube (Carbopack C) | N6200150 |
| TurboMatrix HS Trap Air Monitoring Trap* | M0413628 |